


Iso-Prep Inline Filter System Instructions

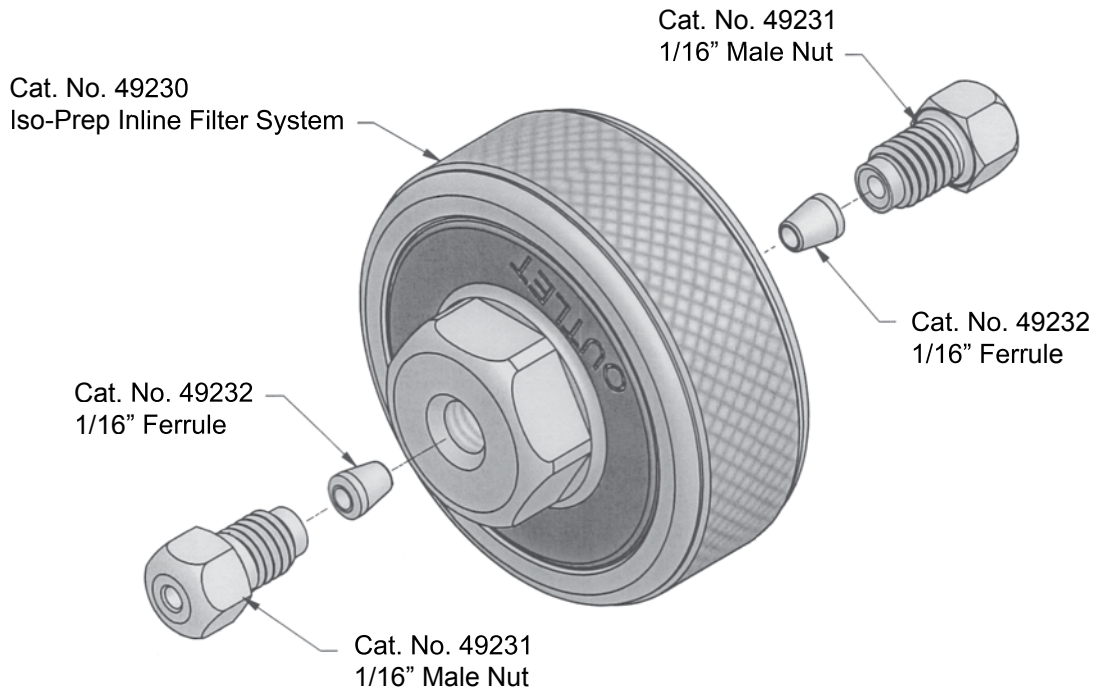

- ▶ Please note: The Iso-Prep Filter Holder has been assembled to a fingertight condition. Using a 9/16" wrench, be sure to torque the Iso-Prep Filter Holder to 40 in/lbs. prior to use.
- ▶ Before use, change existing 1/16" male nuts and ferrules to the supplied male nuts and ferrules to ensure proper connection to the Iso-Prep Filter Holder. **CAUTION!!** Avoid overtightening the 1/16" male nut and ferrule. This may result in damage to the holder.
- ▶ When replacing the filter element inside the Iso-Prep Filter Holder, at re-assembly, re-torque the Iso-Prep Filter Holder between 40 and 60 in/lbs.

Filter replacement: Periodic replacement of the filter element is required to maintain adequate filtration. Filter element lifetime will vary depending on the application. Replacement filter elements (Cat. No. 49234) can be ordered from Analytical Sales and Products.

O-ring replacement: Periodic replacement of the o-ring is required to maintain a proper seal. O-ring lifetime will vary depending on the application. Replacement o-rings (Cat. No. 49233) can be ordered from Analytical Sales and Products.

To order replacement items, please call 973-616-0700.