



Thermal Transfer Deck & Lumidox[®] II LED Array Attachment & Removal Guide

Attaching the Thermal Transfer Deck to a Lumidox® II Solid Base LED Array:

ATTENTION: THIS PROCESS REQUIRES A HIGH LEVEL OF MECHANICAL SKILL TO COMPLETE



STEP 1: Turn the Lumidox II LED array over, and using a 1/16" Allen key, remove the 4 screws securing the bottom plate.



STEP 2: Place the Thermal Transfer Deck on top of the array, and secure it using the 4 longer screws (provided), using the same 1/16" Allen key.



NOTE: Ensure that the screw with the washer is used for the shallow hole near the cable exit location.



STEP 3: Turn the Lumidox II array / Thermal Transfer Deck over. The product is now ready for use.

Removing the Thermal Transfer Deck from a Lumidox® II Solid Base LED Array:

ATTENTION: THIS PROCESS REQUIRES A HIGH LEVEL OF MECHANICAL SKILL TO COMPLETE



STEP 1: Turn the Lumidox II LED array with the Thermal Transfer Deck over, and using a 1/16" Allen key, remove the 4 screws securing them together. Gently lift the TTD away.



STEP 2: Using the same 1/16" Allen key, secure the bottom plate of the Lumidox II with the 4 shorter screws.



NOTE: Ensure that the shortest screw is used for the shallow hole near the cable exit location.



STEP 3: Turn the Lumidox II array over. The product is now ready for use.



179 Rt 206 • Flanders, NJ • 07836
Phone: 973-616-0700 • Fax: 973-616-0133
Email: CustomerService@analytical-sales.com
Website: analytical-sales.com